


# VIDAS® LIS


## Detection and Confirmation of *Listeria* spp

Protocol validated by AOAC International (Official Method of Analysis, Method No. 999.06)  
 \*AOAC RI Performance Tested™ Method Certificate No. 981202

### 1. Detection using VIDAS® LIS

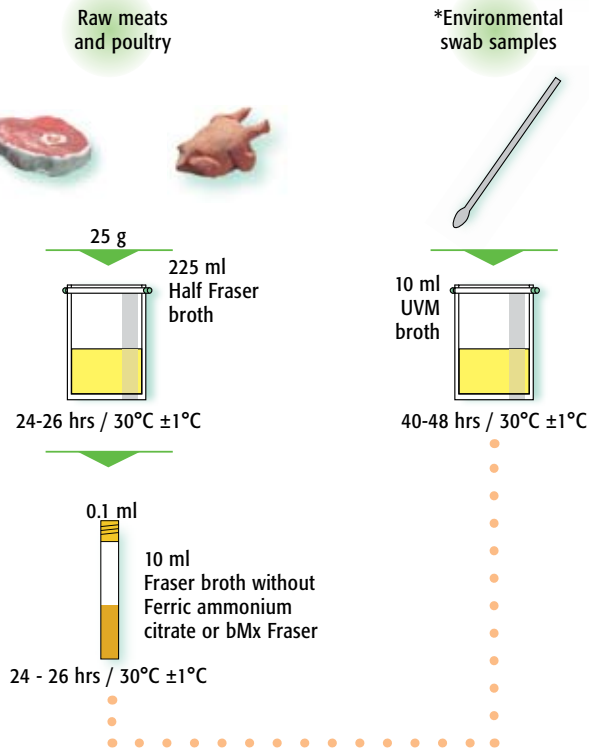

### 2. Confirmation

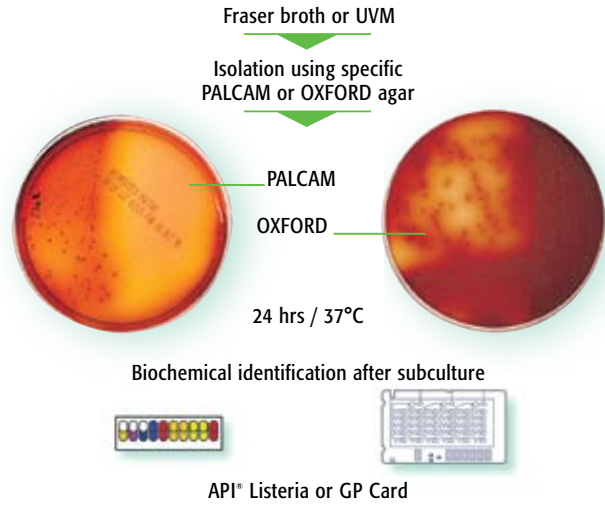

VIDAS® LIS		
7579	Buffered Listeria Enrichment Broth (BLEB)	500 g
T8082	Fraser Broth w/o ferric amm. cit.	10 ml
42046	Fraser Broth	6 x 225 ml
42048	Half Fraser Broth	6 x 225 ml
42072	Fraser Broth	20 x 10 ml
S1210-10	UVM	10 ml
CV 1050	UVM	10 ml
30700	VIDAS LIS	60 tests
M1091	Oxford Agar Modified (MOX)*	10 plates
M1062	LPM Agar	10 plates
7424	LPM Agar	500 g
10300	API Listeria	10 strips
21342	VITEK® 2 GP	20 cards
663141	Listeria Selective Supplement (Oxioid)	

For further details, consult the product package insert.

**bioMérieux, Inc.**  
 595 Anglum Road  
 Hazelwood, MO 63042  
**U.S.A.**  
**Tel: (800) 634 7656**  
**Fax: (800) 657 3053**

**bioMérieux Canada, Inc.**  
 7815 Henri-Bourassa Boulevard West  
 St. Laurent (Québec)  
**Canada H4S 1P7**  
**Tel: (1) 514 336 7321**  
**Fax: (1) 514 807 0015**

[www.biomerieux-usa.com](http://www.biomerieux-usa.com)

