


# VIDAS® LMO 2


## Detection and Confirmation of *Listeria monocytogenes*

Protocol validated by AFNOR & AOAC International (Official Method of Analysis, Method No. 2004.02)  
 \*AOAC approval for all foods products only excluding environments

### 1. Detection using VIDAS® LMO 2

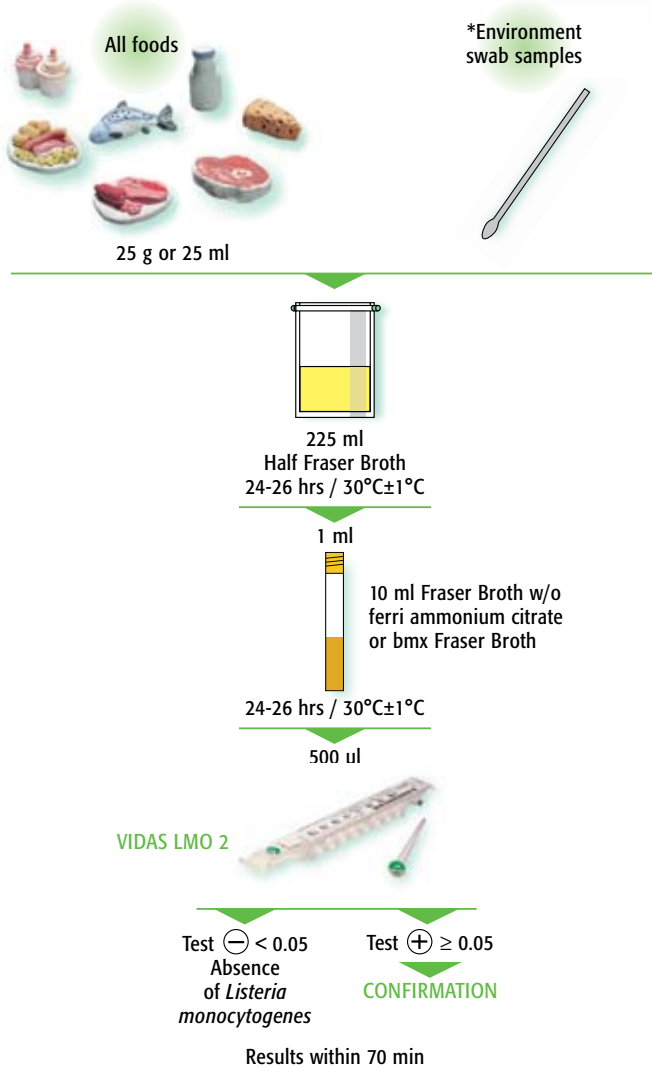

### 2. Confirmation

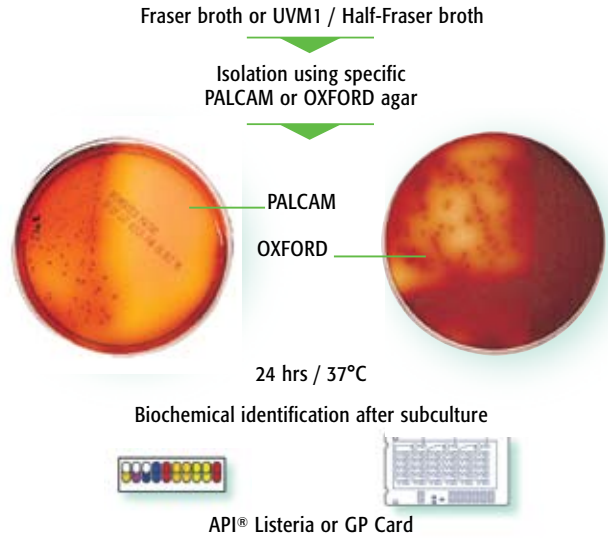

### VIDAS® LMO

42072	Fraser Broth	20 x 10 ml
T8082	Fraser Broth w/o ferric amm. cit.	10 ml
42046	Fraser Broth	6 x 225 ml
S1212-10	UVM	10 ml
CV 1050	UVM	10 ml
42048	Half Fraser Broth	6 x 225 ml
42615	Half Fraser Broth	4 x 225 ml
M1091	Oxford Agar Modified (MOX)*	10 plates
10300	API Listeria	10 strips
30704	VIDAS LMO 2	60 tests
M1062	LPM	10 plates
7424	LPM	500 g
21342	VITEK® 2 GP	20 cards

For further details, consult the product package insert.

**bioMérieux, Inc.**  
 595 Anglum Road  
 Hazelwood, MO 63042  
**U.S.A.**  
**Tel: (800) 634 7656**  
**Fax: (800) 657 3053**

**bioMérieux Canada, Inc.**  
 7815 Henri-Bourassa Boulevard West  
 St. Laurent (Québec)  
**Canada H4S 1P7**  
**Tel: (1) 514 336 7321**  
**Fax: (1) 514 807 0015**

[www.biomérieux-usa.com](http://www.biomérieux-usa.com)

