


# VIDAS® SLM

## Detection and Confirmation of *Salmonella* (TT & SC)

Protocol validated by AOAC International (Official Method of Analysis, Method No. 996.08)



### 1. Detection using VIDAS® SLM

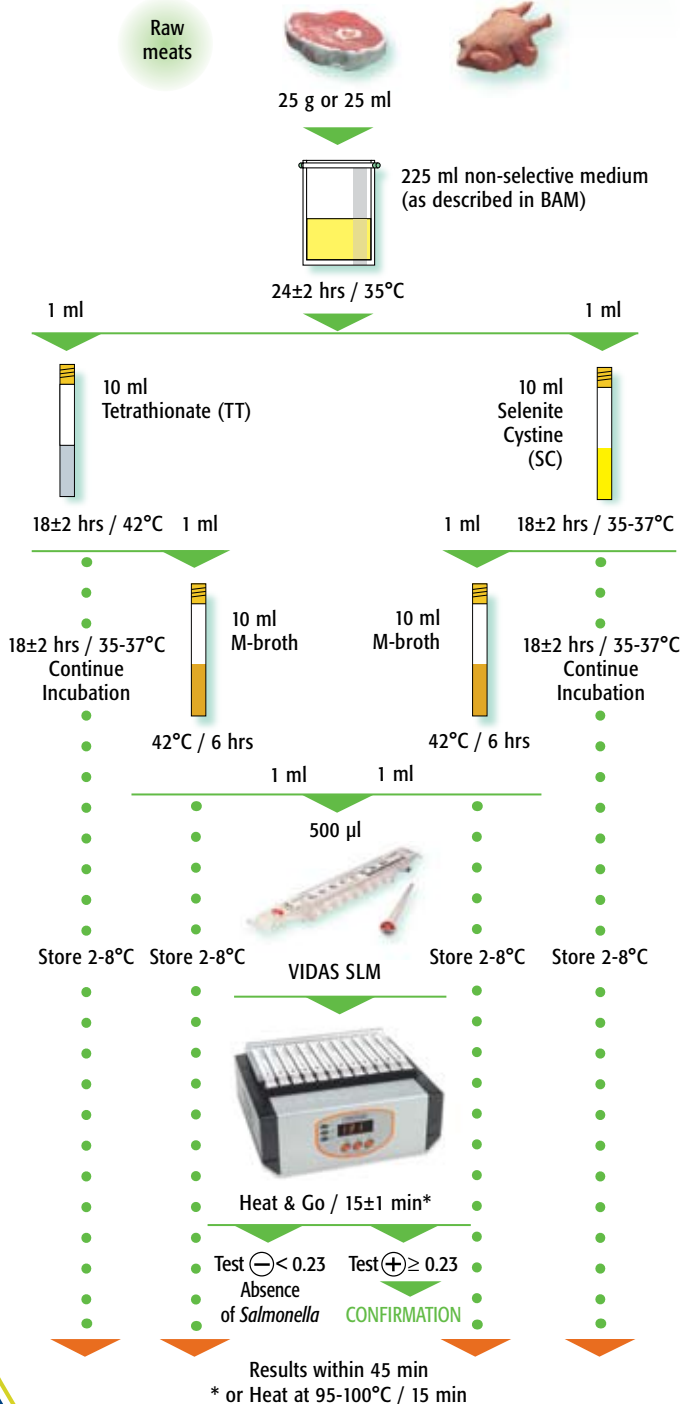

### 2. Confirmation

Streak TT, SC and M broths to Hektoen, XLD and Bismuth Sulfite.  
Incubate plates and confirm suspect colonies

#### Identification



api® System  
For manual  
identification

VITEK® 2 Compact  
For automated  
identification

#### Serological confirmation

#### VIDAS® SLM

T8034	Tetrathionate Broth Base	20 x 10 ml
T7079	Selenite Cystine Broth	20 x 10 ml
T8080	M Broth	20 x 10 ml
30702	VIDAS SLM	60 tests
M1016	Hektoen Enteric Agar	10 plates
M1031	XLD	10 plates
7113	Bismuth Sulfite Agar	500g
43621	ChromID™ Salmonella	20 plates
21341	VITEK 2 GN	20 cards
20100	api 20E	25 strips
32700	rapid ID 32 E	25 strips

For further details, consult the product package insert.

**bioMérieux, Inc.**  
595 Anglum Road  
Hazelwood, MO 63042  
**U.S.A.**  
Tel: (800) 634 7656  
Fax: (800) 657 3053

**bioMérieux Canada, Inc.**  
7815 Henri-Bourassa Boulevard West  
St. Laurent (Québec)  
**Canada** H4S 1P7  
Tel: (1) 514 336 7321  
Fax: (1) 514 807 0015

[www.biomerieux-usa.com](http://www.biomerieux-usa.com)

