

QUANTISWAB™



INNOVATIVE ENVIRONMENTAL MONITORING
MICROBIOLOGY EXPERTISE
UNIQUE & EXCLUSIVE PERFORMANCES




BIOMÉRIEUX
INDUSTRY


QUANTISWAB™




Peel open the 3 wraps to get the sterile QUANTISWAB™ and the tube out

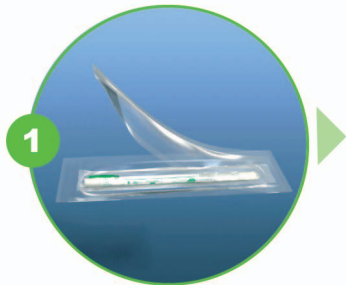

Remove the white cap of the tube

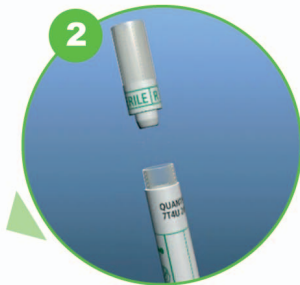

Moisten the QUANTISWAB™ by inserting in the tube and gently squeeze the sponge

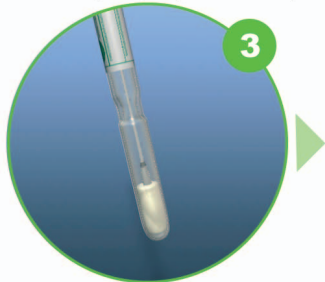

Sample the surface or equipment with QUANTISWAB™



Roll the QUANTISWAB™ directly on an appropriate culture media



Incubate the culture media in the appropriate conditions

